
## MAGUS MCD20/MCD40 MONITOR

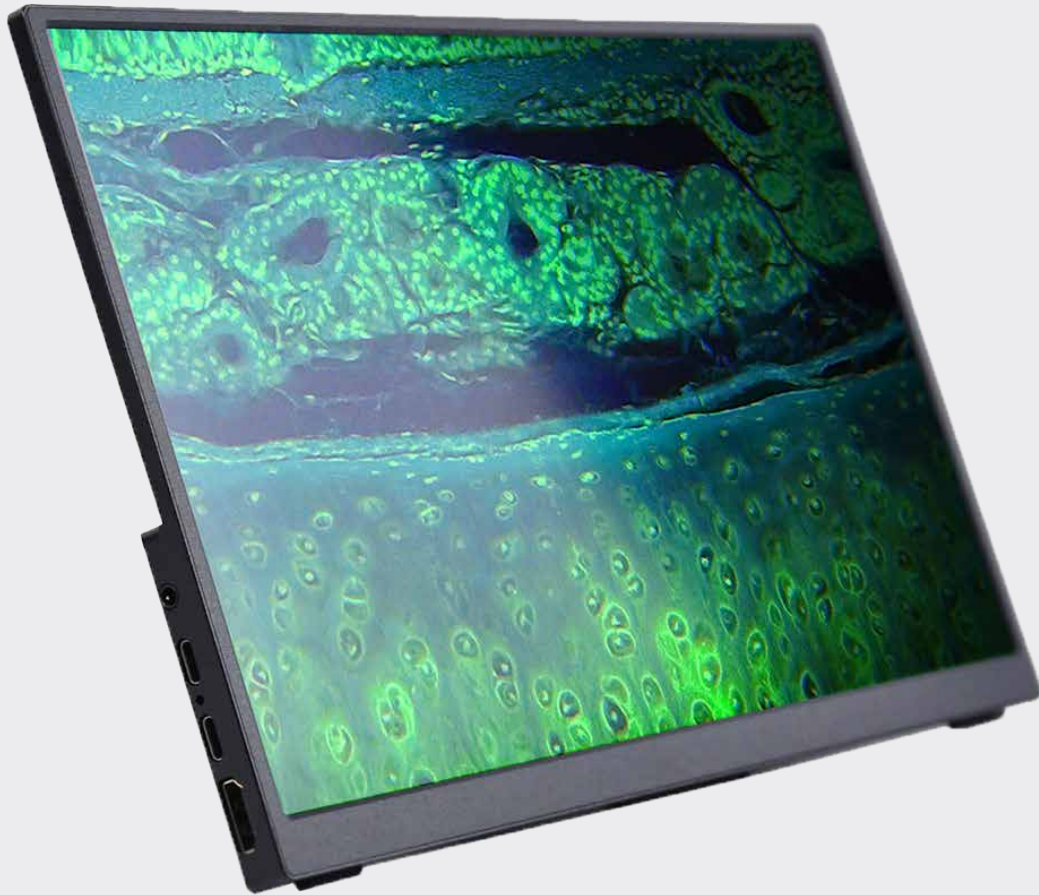# MAGUS

[www.magusmicro.com](http://www.magusmicro.com)

Levenhuk Inc. (USA):  
928 E 124th Ave. Ste D, Tampa, FL 33612, USA,  
+1 813 468-3001,  
[contact\\_us@levenhuk.com](mailto:contact_us@levenhuk.com)

Levenhuk Optics s.r.o. (Europe):  
V Chotejně 700/7, 102 00 Prague 102, Czech Republic,  
+420 737 004-919,  
[sales-info@levenhuk.cz](mailto:sales-info@levenhuk.cz)

Magus® is a registered trademark of Levenhuk, Inc.

© 2006–2024 Levenhuk, Inc. All rights reserved.

## Index

<b>EN</b>	User Manual
<b>BG</b>	Ръководство за потребителя
<b>CZ</b>	Návod k použití
<b>DE</b>	Bedienungsanleitung
<b>ES</b>	Guía del usuario
<b>HU</b>	Használati útmutató
<b>IT</b>	Guida all'utilizzo
<b>PL</b>	Instrukcja obsługi
<b>PT</b>	Manual do utilizador
<b>RU</b>	Инструкция по эксплуатации
<b>TR</b>	Kullanım kılavuzu

# **EN** MAGUS MCD20/MCD40 MONITOR

Before using the monitor, please read this user manual carefully to study the design, operation modes and procedures, operational limitations, and safety precautions.

Due to the continuous improvements in the monitor design, this manual may not reflect minor design changes that do not affect the monitor performance and operation procedures.

## **SAFETY PRECAUTIONS**

Make sure that the input voltage of the monitor matches that of the local power supply. Using the power supply with the wrong input voltage may cause a short circuit or fire.

Do not use a damaged power source.

Do not use a damaged power cord.

Do not insert foreign objects into the slot on the monitor body.

Do not expose the monitor to high temperatures or humidity for a long time.

If water splashes on the monitor, immediately switch the power off, unplug the power cord, and wipe off the water with a dry cloth.

Protect the monitor from shocks, excessive stress from other objects.

Store the monitor away from corrosive environments, household and car heaters, switched-on light bulbs and open flames.

The MAGUS MCD20/MCD40 monitor is designed to create a visualization system of the MAGUS microscope. It connects to a microscope-mounted camera to display a real-time image. The MAGUS MCD20 monitor supports MAGUS Full HD resolution HDMI cameras, while the MCD40 monitor supports 4K resolution cameras.

The monitor is installed on the table or shelf on a folding mount or attached directly to the camera or microscope stand.

The monitor is powered by AC 110–220V, DC 5-12V/1A (Type-C).

# **1 DESCRIPTION OF THE MONITOR**

## **PURPOSE**

The monitor is designed to display a magnified image of the object examined under the microscope. The large resolution provides high image detail and allows you to see the finest details. The IPS matrix delivers a bright image with large viewing angles, which allows you to look at the display even at an angle with no color distortion.

The MAGUS MCD monitors operate in combination with MAGUS CHD cameras to create a complete visualization system for MAGUS microscopes.

The MCD20 reproduces the frame in Full HD resolution and is recommended for MAGUS CHD10, CHD20 and CHD30 cameras.

The MCD40 reproduces the frame in 4K resolution and is recommended for MAGUS CHD40 and CHD50 cameras.

# SPECIFICATIONS

Table 1

	MCD20	MCD40
Type of matrix	IPS	
Screen diagonal, inch	13.3	
Screen resolution, px	1920x1080 (Full HD)	3840x2160 (4K)
Aspect ratio	16:9	
Brightness, cd/m <sup>2</sup>	400	
Number of displayed colors	16.7m	
Contrast	1000:1	
Horizontal/vertical viewing angle, °	178/178	
Size of the visible screen area (WxH), mm	295x165	
Pixel pitch (WxH), mm	0.154x0.154	
Display refresh rate, Hz	60	
Type of matrix backlight	LED	
LED backlight lifetime, h	50000	
Interface	HDMI	
Operating temperature range, °C	-15... 55	
Operating humidity range, %	10... 90	
Power supply	AC 110–220V, DC 5–12V/1A (Type-C)	
Power consumption, W	12 (maximum)	
Package dimensions (WxHxD)	245mm×60mm×355mm	
Weight with package	1.25kg	

The manufacturer reserves the right to make changes to the product range and specifications without prior notice.

# MONITOR KIT

The monitor kit includes the following main components:

- HDMI cable
- AC power adapter
- DC/DC Type-C adapter
- fasteners (Allen wrench and screws)
- User manual and warranty card.

See Section 5 of the User manual for a full kit contents.

The general view of the monitor is given in Fig. 1.

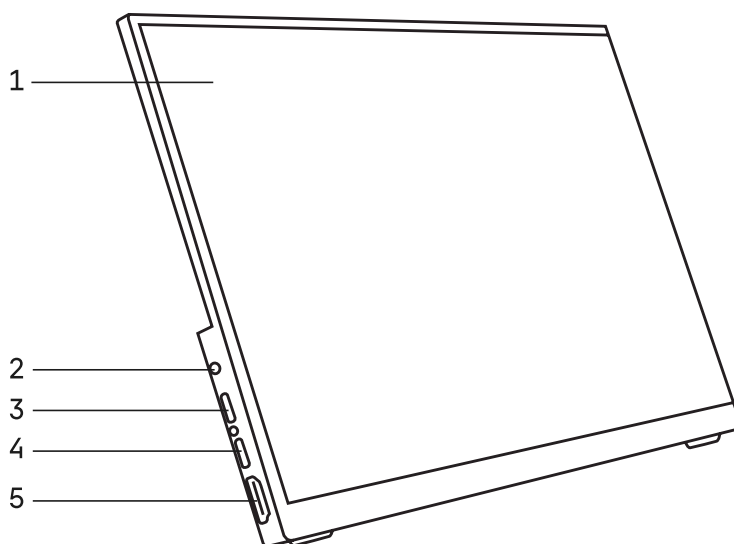
Fig. 1. MCD20/MCD40 Monitor

- |                     |                     |
|---------------------|---------------------|
| 1 – monitor         | 4 – USB Type-C port |
| 2 – audio output    | 5 – HDMI port.      |
| 3 – USB Type-C port |                     |

## 2 UNPACKING THE MONITOR

Unpack the monitor.

Check the scope of delivery using Section 5 of the User manual.

Install the monitor on the table or shelf on a folding mount, or attach it directly to the camera or microscope stand.

Make sure that the monitor is securely and safely installed.

Check and sort the supplied accessories and tools in the correct order. Keep them in proper order to avoid confusion.

### 3 USING THE MONITOR

To display the image on the monitor screen:

Attach the monitor 4 to the camera 1 using the fasteners from the kit.

Connect the monitor to the camera using the HDMI cable 3.

Connect the monitor and camera to AC power using the DC/DC Type-C adapter and power adapter (supplied). If the camera and monitor are remote from each other, each device is powered separately using the power adapter supplied with both the camera and monitor.

Switch on and adjust the camera as described in the manual supplied with the camera.

Switch on the monitor by pressing the bottom button (not shown in Figure 2) on the side panel as indicated by the arrow.

If the image on the screen is blurred, rotate the fine focusing knob to get an accurate image.

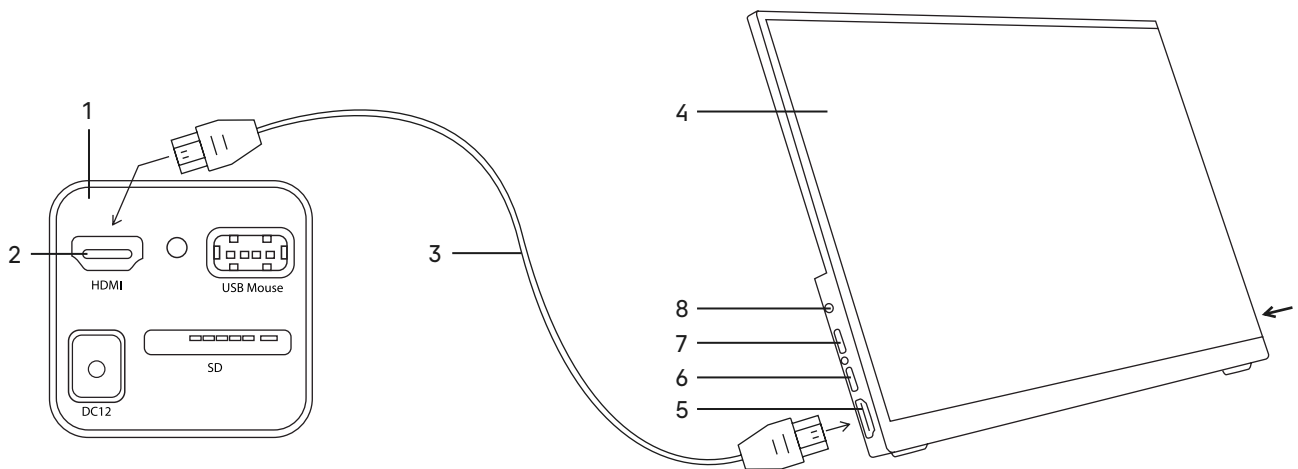
Fig. 2. Using the monitor

- 1 – camera
- 2 – HDMI port
- 3 – HDMI cable

- 4 – monitor
- 5 – HDMI port
- 6 – USB Type-C port

- 7 – USB Type-C port
- 8 – audio output

## 4 TROUBLESHOOTING

Potential problems and remedies are given in Table 2.

Table 2

Problem	Cause	Remedy
There is no image on the monitor screen	The power supply is faulty	Replace the power supply
	The power cord is faulty	Replace the power cord
	The power connector has poor contact	Have the connector repaired by a qualified electronics technician
	The camera is faulty	Have the camera repaired by a qualified electronics technician
The image does not remain sharp during observation	The wrong camera has been selected	Select the appropriate camera
	The focus is not adjusted	Adjust the focus using the focusing knob
The specimen image is not in the center of the field of view	The camera position is not adjusted	Adjust the camera position

## 5 SCOPE OF DELIVERY

Full scope of delivery (Table 3)

Component	Pcs	Note
Monitor	1	
HDMI cable	1	
AC power adapter	1	
DC/DC Type-C adapter	1	
Fasteners (Allen wrench and screws)	1	
User manual and warranty card	1	

## 6 MAINTENANCE OF THE MONITOR

1. Once you have finished using the monitor, switch off the power supply. When not using the monitor for a long time, unplug the power cord from the power outlet.

2. Store the monitor in a dry, ventilated and clean place, with no exposure to acids, alkalis, or steam.

3. The monitor should be kept clean. Use a soft brush to remove dust and small particles. If there are liquid drops on the screen, remove them with a dry cloth or soft tissue. Use special alcohol wipes to remove heavy soiling.

**Always switch off the monitor before cleaning. Do not use aggressive agents to clean difficult stains, as this may cause damage to the device.**

4. Periodic inspection: the monitor should be regularly inspected and serviced to maintain its performance

## 7 MAGUS WARRANTY

MAGUS provides a **5-year international warranty** from date of purchase (valid for the entire life of the instrument). The Levenhuk company warrants the product to be free from defects in materials and workmanship. The Seller warrants that the MAGUS product you have purchased meets specification requirements, provided that the Buyer complies with terms and conditions of transport, storage, and operation of the product. The warranty period for accessories is **6 (six) months** from the date of purchase.

For more information on warranty terms and conditions, see [www.magusmicro.com](http://www.magusmicro.com)

For warranty service, please contact your nearest Levenhuk representative office.

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